

MC6000 Replacement Syringes

| | | | | | | | |
|---------------|------------|-------------|-------------|-------------|-------|--------|-------|
| Cat. # | 10080 | 10081 | 10082 | 10083 | 10084 | 10085 | 10086 |
| Volume | 50 μ L | 100 μ L | 250 μ L | 500 μ L | 1 mL | 2.5 mL | 5 mL |

***Materials:** PTFE seal, Borosilicate Glass

Fitting Interface: 1/4 -28

Stroke: 3 cm

*UHMWPE Syringes are available upon request.



SYRINGE REPLACEMENT

To replace the syringe, prime all fluid from the syringe, and follow the steps below:

1. To replace the syringe, the syringe plunger holder must be in the lowered position. This can be done by sending the A3000R command.
2. Turn the plunger lock screw counterclockwise 3 turns.
3. Turn the syringe counterclockwise and remove from the valve.

To install the syringe follow the steps below:

4. Screw the threaded portion of the syringe clockwise into the valve. Align barrel plunger to center of hole in the plunger holder.
5. Fully seat the plunger into the syringe plunger holder.
6. Securely tighten (clockwise) the plunger lock screw.
7. Initialize the pump.

